



Cryogenic Workwear Ltd
Protective equipment for cryogenic operatives

Tel: +44 (0)1908 221755
www.cryogenic-gloves.co.uk

Face Shield

LENS MARKING - 2C-1.2 1 AT 8 9 CE

FRAME MARKING - EN166 3 8 9 AT CE



High quality easy to fit face shield, with height and width adjustable headgear and headband.

The injection moulded translucent polycarbonate shield is shaped to hug the side of your face, which offers good protection against liquid droplets and vapour, and is the strongest material for impact resistance, but is also lightweight.

This face shield and visor gives a 180° field of unrestricted vision, whilst providing facial protection.

A flip up shield is incorporated to enable the user to lift up, when their work has been completed.

Visors should be inspected before each use to ensure that they are not damaged. They can be cleaned using warm soap and water, the use of harsh abrasives is not recommended.

In use, ensure that the face shield fits comfortably and securely on the head. Protection will only be offered if the face shield is fitted correctly.

Ensure that the visor is in place before use.

The face shield should be stored out of direct sunlight, away from chemicals and not exposed to extremes of temperature.